

 **BD Rhapsody™ System**  
Single-Cell Labeling with  
BD® AbSeq Ab-Oligos (1 plex to 40 plex)  
Protocol

## Copyrights

No part of this publication may be reproduced, transmitted, transcribed, stored in retrieval systems, or translated into any language or computer language, in any form or by any means: electronic, mechanical, magnetic, optical, chemical, manual, or otherwise, without prior written permission from BD.

The information in this guide is subject to change without notice. BD reserves the right to change its products and services at any time. Although this guide has been prepared with every precaution to ensure accuracy, BD assumes no liability for any errors or omissions, nor for any damages resulting from the application or use of this information. BD welcomes customer input on corrections and suggestions for improvement.

## Trademarks

BD, the BD Logo, BD Rhapsody and Fc Block are trademarks of Becton, Dickinson and Company or its affiliates. All other trademarks are the property of their respective owners. © 2022 BD. All rights reserved.

For US patents that may apply, see [bd.com/patents](https://bd.com/patents).

## Regulatory information

For Research Use Only. Not for use in diagnostic or therapeutic procedures.

## History

Revision	Date	Change made
Doc ID: 214394 Rev. 1.0	2018-07	Initial release.
23-24262(01)	2022-11	Added BD Rhapsody™ Enhanced Cell Capture Beads v2.0. Removed part numbers.

# Contents

---

<b>Introduction</b> .....	<b>4</b>
<b>Required materials</b> .....	<b>4</b>
<b>Suggested materials</b> .....	<b>4</b>
<b>Before you begin</b> .....	<b>4</b>
Safety information .....	5
<b>Preparing 2X BD® AbSeq antibody-oligo labeling MasterMix</b> .....	<b>5</b>
<b>Labeling cells with BD® AbSeq Ab-Oligos</b> .....	<b>6</b>
<b>Washing labeled cells</b> .....	<b>7</b>
<b>Troubleshooting</b> .....	<b>8</b>

## Introduction

This protocol describes use of BD® AbSeq Ab-Oligos (antibody-oligonucleotides) for antigen-expression profiling with BD Rhapsody™ single-cell capture and downstream library preparation. Each BD® AbSeq Ab-Oligo is an oligonucleotide-conjugated antibody that contains an antibody-specific barcode and poly(A) tail for bead capture, PCR amplification, and library generation. The protocol supports BD® AbSeq Ab-Oligo labeling of 20,000 to 1 million cells. Up to 40 antibodies can be pooled together per staining reaction. For greater than 40 antibodies, refer to *BD Rhapsody™ System Single-Cell Labeling with BD® AbSeq Ab-Oligos (41 plex to 100 plex) Single-Cell Protocol*.

## Required materials

Material	Supplier	Catalog no.
20,000-1 million cells	–	–
BD® Stain Buffer (FBS)	BD Biosciences	554656
BD® AbSeq Ab-Oligos <sup>a</sup>	BD Biosciences	Various
BD Rhapsody™ Enhanced Cartridge Reagent Kit	BD Biosciences	664887
Latch Rack for 500-µL tubes	Thermo Fisher Scientific	4900 or 4890
Falcon® tubes, 5-mL round-bottom, polystyrene test tubes <sup>b</sup>	Corning	352054
a. Avoid storing BD® AbSeq Ab-Oligos or Sample Tags under freezing conditions.		
b. Use only the tubes specified in the protocol. Use of other tubes might lead to increased cell loss.		

For a complete list of materials, see appropriate instrument user guide.

## Suggested materials

Material	Supplier	Catalog no.
Human BD Fc Block™	BD Biosciences	564220
Or,		
Mouse BD Fc Block™	BD Biosciences	553142
8-Channel Screw Cap Tube Capper	Thermo Fisher Scientific	4105MAT
Multi-channel pipette	Major supplier	–

## Before you begin

- Use low retention filtered pipette tips.
- Prepare a single-cell suspension. See *Preparing Single-Cell Suspensions Protocol*.
- If your biological sample contains red blood cell contamination, red blood cell lysis is required. See *Preparing Single-Cell Suspensions Protocol*.

## Safety information

For safety information, see the *BD Rhapsody™ Single-Cell Analysis Instrument User Guide* or the *BD Rhapsody™ Express Single-Cell Analysis System Instrument User Guide*.

## Preparing 2X BD® AbSeq antibody-oligo labeling MasterMix

We recommend:

- Creating freshly pooled antibodies before each experiment.
  - Creating pools with 30% overage to ensure adequate volumes for labeling. The reagents are viscous and form bubbles easily.
  - For high-plex panels, using an 8-Channel Screw Cap Tube Capper and multi-channel pipette to pipet BD® AbSeq Ab-Oligos into 8-tube strips. Centrifuge tube strip and pool BD® AbSeq Ab-Oligos into a 1.5-mL LoBind tube.
- 1 Place all BD® AbSeq Ab-Oligos to be pooled into a Latch Rack for 500 µL tubes. Arrange the tubes so that they can be easily uncapped and re-capped with an 8-Channel Screw Cap Tube Capper and aliquoted with a multi-channel pipette.
  - 2 Centrifuge BD® AbSeq Ab-Oligos in the Latch Rack in a tabletop centrifuge with a plate adapter tubes at  $400 \times g$  for 30 seconds and place on ice.
  - 3 In pre-amplification workspace, pipet reagents into a new 1.5-mL LoBind Tube on ice:

### 2X BD® AbSeq labeling MasterMix

Component	1 sample (µL)	1 sample + 30% overage (µL)	2 samples + 30% overage (µL)
Per BD® AbSeq Ab-Oligo	2.0	2.6	5.2
BD Stain Buffer (FBS) ( <i>N</i> = no. antibodies)	100 – (2.0 × <i>N</i> )	130 – (2.6 × <i>N</i> )	260 – (5.2 × <i>N</i> )
<b>Total</b>	<b>100.0</b>	<b>130.0</b>	<b>260.0</b>

### Examples

Component	1 sample (µL)	1 sample + 30% overage (µL)	2 samples + 30% overage (µL)
<b>10-plex BD® AbSeq labeling</b>			
Per BD® AbSeq Ab-Oligo	2.0 (20.0 total)	2.6 (26.0 total)	5.2 (52.0 total)
BD Stain Buffer (FBS)	80.0	104.0	208.0
<b>20-plex BD® AbSeq labeling</b>			
Per BD® AbSeq Ab-Oligo	2.0 (40.0 total)	2.6 (52.0 total)	5.2 (104.0 total)
BD Stain Buffer (FBS)	60.0	78.0	156.0
<b>40-plex BD® AbSeq labeling</b>			
Per BD® AbSeq Ab-Oligo	2.0 (80.0 total)	2.6 (104.0 total)	5.2 (208.0 total)
BD Stain Buffer (FBS)	20.0	26.0	52.0

- 4 Pipet-mix the 2X BD® AbSeq labeling MasterMix, and place back on ice.

## Labeling cells with BD® AbSeq Ab-Oligos

- 1 Centrifuge cells at  $400 \times g$  for 5 minutes, or at an appropriate speed to pellet the cells.
- 2 (Optional) For samples containing myeloid and B lymphocytes, we recommend blocking non-specific Fc Receptor-mediated false-positive signal with Human BD Fc Block™ or Mouse BD Fc Block™, as appropriate.

To perform blocking:

- a Pipet reagents into a new 1.5-mL LoBind tube on ice:

### Fc Block™ MasterMix

Component	For 1 sample (μL) <sup>a</sup>	For 1 sample + 20% overage (μL)
BD Stain Buffer (FBS)	95.0	114.0
Human BD Fc Block™ or Mouse BD Fc Block™	5.0	6.0
<b>Total</b>	<b>100.0</b>	<b>120.0</b>

a. Sufficient for  $\leq 1 \times 10^6$  cells. To block more cells, adjust volume.

- b Pipet-mix Fc Block™ MasterMix and briefly centrifuge. Place on ice.
  - c Discard supernatant from cells without disturbing pellet.
  - d Resuspend cells in 110 μL Fc Block™ MasterMix.
  - e Incubate cells at room temperature (15–25 °C) for 10 minutes.
  - f After Fc Block™, proceed to **step 4**.
- 3 Discard supernatant from cells without disturbing pellet, and resuspend each sample in 110 μL BD® Stain Buffer. Pipet-mix.
  - 4 In a new 5-mL polystyrene Falcon® tube, combine 100 μL of cell suspension and 100 μL 2X BD® AbSeq labeling MasterMix. Pipet-mix
  - 5 Incubate on ice for 30–60 minutes.

## Washing labeled cells

---

**Note:** Sufficient post-labeling washing is important for reducing noise that comes from residual unbound antibodies being captured onto 3' capture beads during single-cell capture. However, some cell loss occurs with each additional wash. Users can choose to perform more or fewer washes depending on the abundance of their sample.

- 1 Add 2 mL BD® Stain Buffer to labeled cells and pipet-mix.
- 2 Centrifuge each tube at  $400 \times g$  for 5 minutes, or at an appropriate speed to pellet the cells.
- 3 Uncap each tube and invert to decant supernatant into biohazardous waste. Keep the tube inverted and gently blot on a lint-free wiper to remove residual supernatant from tube rim.
- 4 Add 2 mL BD® Stain Buffer to each tube, and resuspend by pipet-mixing.
- 5 Centrifuge at  $400 \times g$  for 5 minutes, or at an appropriate speed to pellet the cells.
- 6 Uncap each tube, and invert to decant supernatant into biohazardous waste. Keep the tube inverted and gently blot on a lint-free wiper to remove residual supernatant from tube rim.
- 7 (Optional) Repeat **steps 4–6** once more for a total of 3 washes.
- 8 Resuspend the pellet in 620  $\mu$ L cold Sample Buffer from the BD Rhapsody™ Enhanced Cartridge Reagent Kit. Perform viability staining and count cell using the appropriate single-cell capture and cDNA synthesis protocol.  
  
**Note:** For low-abundance samples (<20,000), resuspend the cells in 200  $\mu$ L of cold BD® Sample Buffer. For other 3' single-cell capture platforms, resuspend in recommended buffer and volume according to manufacturer.
- 9 Place tube on ice, and proceed to single-cell capture. See the Single-Cell Analysis Workflow with BD Rhapsody™ Systems to find the appropriate protocol to follow.

## Troubleshooting

Observation	Possible causes	Recommended solutions
Do not have the recommended buffer for labeling with BD <sup>®</sup> AbSeq Ab-Oligos.	Various.	Labeling with BD <sup>®</sup> AbSeq Ab-Oligos is optimal in BD <sup>®</sup> Stain Buffer (FBS). Label BD <sup>®</sup> AbSeq Ab-Oligos in BD <sup>®</sup> Stain Buffer (FBS).
Cells require labeling with BD <sup>®</sup> AbSeq Ab-Oligos at a different temperature.	Physiological requirement	Use protocols for BD <sup>®</sup> AbSeq Ab-Oligo labeling that have been optimized for the specific sample type.
Accidentally resuspended cells in BD <sup>®</sup> Stain Buffer (FBS) rather than Sample Buffer before cell counts.	Various.	We recommend centrifuging the samples and resuspending the cells in Sample Buffer after labeling with BD <sup>®</sup> AbSeq Ab-Oligos. This ensures optimal performance of cell loading in the BD Rhapsody™ Cartridge.
Cell loss.	Wrong tube used in washes.	Use Falcon <sup>®</sup> polystyrene flow tubes and centrifuge cells using a benchtop centrifuge with swing bucket rotor. This centrifugation method reduces cell loss.
Cell loss after sorting.	Various.	<ul style="list-style-type: none"> <li>Sort more cells than needed for cartridge loading.</li> <li>Sort cells into 5 mL polystyrene Falcon<sup>®</sup> tube. Use the same 5 mL polystyrene Falcon<sup>®</sup> tube that was used for sorting for cell labeling by following these steps:               <ol style="list-style-type: none"> <li>Create a 1X AbSeq labeling MasterMix by adding 100 <math>\mu</math>L BD<sup>®</sup> Stain Buffer per 100 <math>\mu</math>L 2X BD<sup>®</sup> AbSeq labeling MasterMix.</li> <li>Pipet-mix, and place on ice.</li> <li>Sort cells into a 5 mL polystyrene Falcon<sup>®</sup> tube.</li> <li>Centrifuge the sorted cell suspension at 400 <math>\times</math> g for 5 minutes.</li> <li>Uncap the tube and invert to decant supernatant into biohazardous waste.</li> <li>Keep the tube inverted and gently blot on a lint-free wiper to remove residual supernatant from tube rim.</li> <li>Resuspend cell pellet with the 1X BD<sup>®</sup> AbSeq labeling MasterMix (<b>step 1</b>), and proceed with cell labeling.</li> </ol> </li> </ul>



**Becton, Dickinson and Company**  
**BD Biosciences**  
2350 Qume Drive  
San Jose, California 95131 USA

[bdbiosciences.com](http://bdbiosciences.com)  
[scomix@bdscomix.bd.com](mailto:scomix@bdscomix.bd.com)